


Mounting instruction for the fill hose to the pump

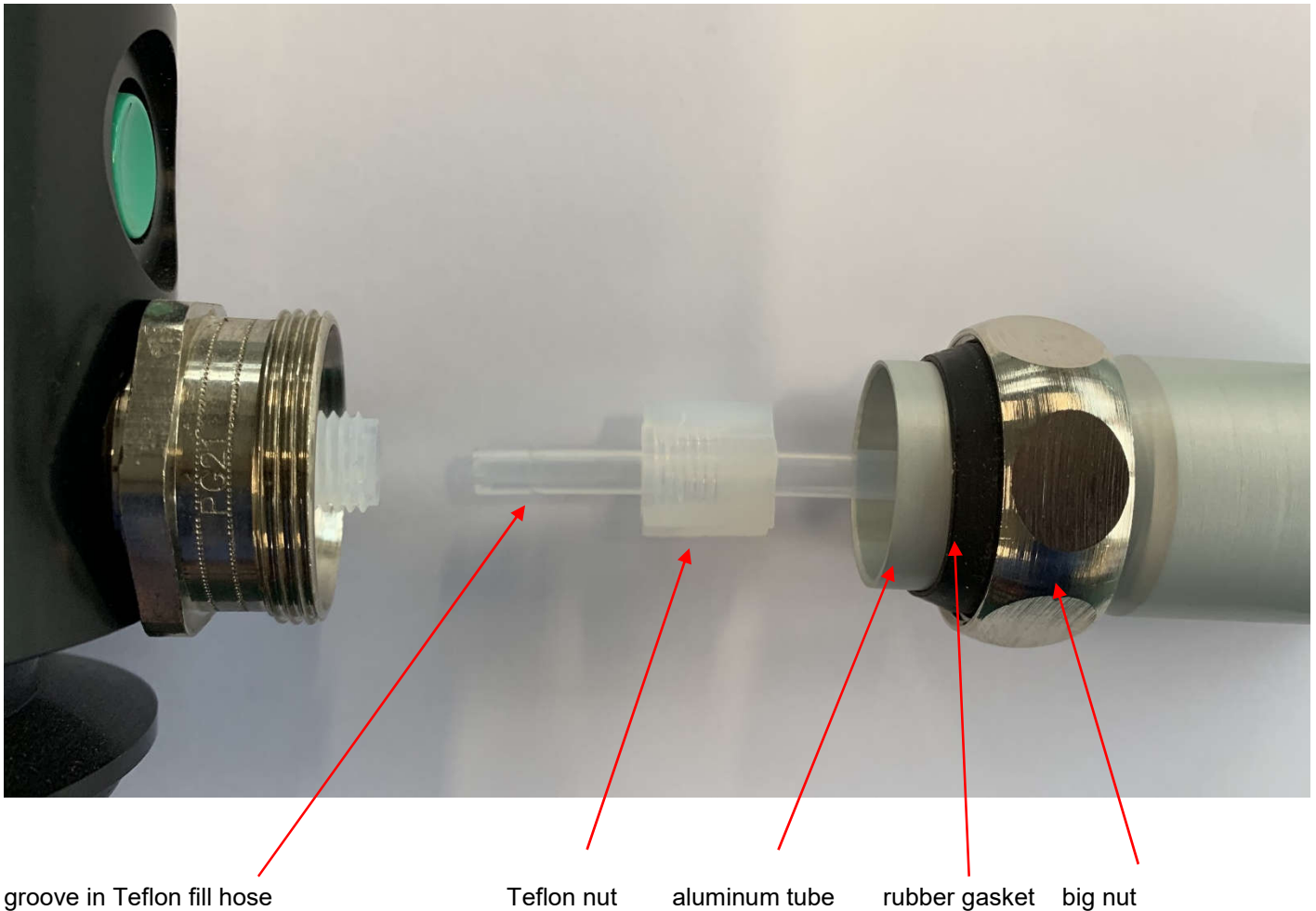

- Loosen the big nut on the outlet, and put it on the metal protection tube of the fill hose. Also slide the black rubber gasket as far over the aluminum tube as possible. Loosen the Teflon (plastic) nut 2 turns. Push the Teflon (plastic) fill hose in the nut. Tighten the Teflon nut 2 turns, and **gently pull on the fill hose until the groove clicks in the nut**. This is easy to feel. Now tighten the nut by hand as tight as possible. Then use a spanner to give the nut one full turn. DO NOT TIGHTEN IT TOO MUCH. It is plastic, and can break if tightened too much.
- Move the metal protection tube into the pump's outlet. Tighten the big nut so that the protection tube is fixed. Tighten by hand only.

For a clear instruction how to Leak Free install the Cryo hose, please check our Norhof YouTube instruction Video at:
https://youtu.be/xy8xilgeiJY?si=eplJ_bDJgx2Qbslk

